

# dataLink™

Traceability system for easySpiral® platers  
and Scan® automatic colony counters

## 100% traceability

from plating to colony counting



interscience



easy**Spiral** Dilute®  
serial dilutor and automatic plater



**dataLink™** is a full traceability system for the microbiological analysis of a sample, from plating to colony counting with **easySpiral®** automatic platers and **Scan®** automatic colony counters.

All plating data (volume, dilution, time, mode ...) is stored on a Datamatrix or QR Code label placed on the Petri dish. No database is needed.

After incubation, the Datamatrix label is scanned and read by the **Scan®** software and retrieves all the sample information, from the plating to the colony counter settings.

Simply press "COUNT". There is no more time-consuming data double input and potential errors. Export the results, pictures and comments to LIMS\*, Excel™, PDF, JPEG.

### ▶ **Secured analyses**

- Full traceability
- Error-free settings

### ▶ **Time-saving**

- No manual data input
- Automatic settings of the Scan® colony counter

### ▶ **User-friendly**

- Stand-alone system
- No database or network use

# STEP 1

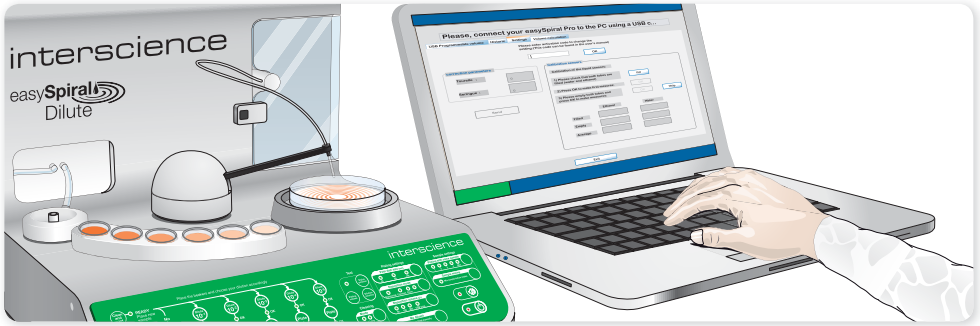


Plate with easySpiral® Pro or Dilute. easySpiral® software collects the plating data.

# STEP 2



Print the label with Datamatrix code. Stick the label on the plated Petri dish and place in the incubator.

■■■ Incubation 24-72 h

# STEP 3



Once the colonies have grown, scan the Datamatrix code. The **Scan®** colony counter automatically adjusts its settings thanks to the Datamatrix label's data. Click on "COUNT". Export the data.

\* Please check LIMS compatibility

# 5 label models:




## Bar-code 39



## Bar-code 128



## Datamatrix

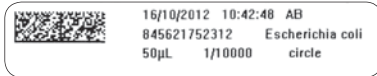

## QR Code



## Number



Label dimensions: 45 x 10 mm

# dataLink™ pack:



Ref. 410 100



Your local distributor

## interscience INTERNATIONAL

30, ch. Bois Arpents - 78860 St Nom - FRANCE  
T: +33 (0)1 34 62 62 61 - F: +33 (0)1 34 62 43 03  
E mail: info@interscience.com

## interscience USA & CANADA

32 Cummings Park - Woburn, MA 01801 - USA  
T: +1 781 937 0007 - F: +1 781 937 0017  
E mail: sales.usa@interscience.com

## interscience CHINA

上海徐汇区吴兴路277号锦都大厦718室-200030  
电话: +86(0)21-64739390 - 手机: +86 189 3097 0733  
邮址: sales.china@interscience.com

## interscience SOUTH-EAST ASIA

541 Orchard Rd -09-01 Liat Towers - SINGAPORE 238881  
T: +65 6933 1389 - M: +65 8549 1217  
E mail: sales.asia@interscience.com